

# NEUROMICS



## Cortical Neurons Kit

---

**Catalog #:** PC35117

**Product Format:** T25 Flask

---

Enclosed is a T25 culture flask containing live adherent primary cortical neurons obtained from Sprague Dawley rats and a 12 ml tube of NbActiv4® culture medium (Neurobasal®/B27®/GlutaMAX™). Live adherent neurons are stored in ~65ml of our shipping medium. For cell survival, an immediate media exchange is required.

Preparations:

1. Remove the T25 culture flask and 12 ml tube of NbActiv4® medium from the shipping container
2. Allow both the T25 culture flask and the 12 ml tube of NbActiv4® medium to equilibrate to room temperature

Media Exchange (Room Temperature in a Sterile Hood)

1. Remove the parafilm seal
2. Sterilize the T25 culture flask lid with 70% EtOH and allow to air dry
3. For each T25 culture flask (one flask at a time)
4. Aspirate off the shipping media
5. Immediately add 5.0 mL of NbActiv4® medium to the T25 culture flask
6. Repeat for all remaining T25 flasks
7. Incubate the T25 culture flask(s) at 37°C, 5% CO<sub>2</sub>, 9% O<sub>2</sub>, 95% humidity (or ambient O<sub>2</sub>)
8. Upon receipt, neurons display axons and dendrites; synapses and action potentials begin at 7 days
9. Change ½ of the medium with fresh, 37°C, CO<sub>2</sub> equilibrated NbActiv4® every 3-4 days.

Viability Assay:

1. Rinse twice with 37°C HBSS (0.2 ml/cm<sup>2</sup> of substrate).
2. Prepare dye mix from an acetone stock of 15 mg/ml fluorescein diacetate and an aqueous stock of 4.6 mg/ml propidium iodide, dilute 15 µl of each into 1.5 ml HBSS (1:100 dilution).
3. Add 20 µl of dye mix from step 2 to every 0.2 ml of HBSS added in step 1 (1:10 dilution)
4. Afer ~1 minute count live cells using blue excitation appropriate for fluorescein fluorescence (green cells). Count dead cells with green excitation for propidium iodide fluorescence (small red nuclei)
5. Viability = (green cells/unit area)/(total cells plated/unit area) or Survival = green cells/(green + red cells)

Neurobasal®/B27®/GlutaMAX™ are trademarks of Life Technologies.

### FOR RESEARCH USE ONLY

NEUROMICS REAGENTS ARE FOR IN VITRO AND CERTAIN NON-HUMAN IN VIVO EXPERIMENTAL USE ONLY AND NOT INTENDED FOR USE IN ANY HUMAN CLINICAL INVESTIGATION, DIAGNOSIS, PROGNOSIS, OR TREATMENT. THE ABOVE ANALYSES ARE MERELY TYPICAL GUIDES. THEY ARE NOT TO BE CONSTRUED AS BEING SPECIFICATIONS. ALL OF THE ABOVE INFORMATION IS, TO THE BEST OF OUR KNOWLEDGE, TRUE AND ACCURATE. HOWEVER, SINCE THE CONDITIONS OF USE ARE BEYOND OUR CONTROL, ALL RECOMMENDATIONS OR SUGGESTIONS ARE MADE WITHOUT GUARANTEE, EXPRESS OR IMPLIED, ON OUR PART. WE DISCLAIM ALL LIABILITY IN CONNECTION WITH THE USE OF THE INFORMATION CONTAINED HEREIN OR OTHERWISE, AND ALL SUCH RISKS ARE ASSUMED BY THE USER. WE FURTHER EXPRESSLY DISCLAIM ALL WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. v1-09809

[www.neuromics.com](http://www.neuromics.com)

Neuromics Antibodies • 5325 West 74<sup>th</sup> Street, Suite 8 • Edina, MN 55439  
Phone 866-350-1500 • fax 612-677-3976 • email: [pshuster@neuromics.com](mailto:pshuster@neuromics.com)