

Data Sheet

Gravity flow Strep-Tactin®XT 4Flow® **High Capacity column**

Cat. No.: 2-5031-001, 2-5031-005, 2-5032-001,

2-5033-001, 2-5034-001

Lot No.:

IBA Headquarters IBA GmbH

Rudolf-Wissell-Str. 28 37079 Goettingen Germany Tel. +49 (0) 551-5 06 72-0 Fax +49 (0) 551-5 06 72-181

IBA US contact information

Fax 1-888-531-6813

Version:

E-mail: info@iba-lifesciences.com https://www.iba-lifesciences.com

1.0 Revision Date: 02.03.2020

Description	Ready-to-use column with Strep-Tactin®XT high capacitiy resin for the purification of Strep-tag®II and Twin-Strep-tag® fusion proteins. Strep-Tactin®XT is a streptavidin variant with optimized binding properties for Strep-tag® fusion proteins.
Support	4% agarose, crosslinked
Form	Pre-packed in buffer, pH 8.0: 100 mM Tris/HCl pH 8.0, 1 mM EDTA, 150 mM NaCl, 0.02% sodium azide
Dynamic Binding Capacity	16 mg protein/ml resin. Dynamic binding capacity was determined with 1 mg/ml mCherry-Twin-Strep-tag® (30 kDa) at a flow rate of 0.5 ml/min. Please note: Binding capacity is protein dependent.
Stability	6 months after shipping
Storage	recommended: 2–8 °C
Shipping	room temperature
Hazards	Product is not classified as hazardous according to (EC) No 1272/2008 [CLP]. A Material Safety Data Sheet is provided.

Elution	Strep-Tactin®XT Elution Buffer with Biotin (10 x Buffer BXT, Cat. No. 2-1042-025); 1 x Buffer BXT: 100 mM Tris-HCl pH 8.0, 150 mM NaCl, 1 mM EDTA, 50 mM biotin
Regeneration	Regenerate the column by using Strep-Tactin®XT Regeneration Buffer (3 M MgCl ₂ , Cat. No. 2-1045-250)

Voss, S. & Skerra, A. (1997) Mutagenesis of a flexible loop in streptavidin leads to higher affinity for the Strep-tag II peptide and improved performance in recombinant protein purification. Protein Eng. 10, 975-982.



Please download all up-to-date manuals, protocols and other material from https://www.iba-lifesciences.com.

For research use only

This product is based on Strep-Tactin®XT technology covered by intellectual property (IP) rights and on completion of the sale IBA grants respective Limited Use Label Licenses to purchaser. IP rights and Limited Use Label Licenses for said technology are further described and identified at http://www.iba-lifesciences.com/patents.html or upon inquiry at info@iba-lifesciences.com or at IBA GmbH, Rudolf-Wissell-Str. 28, 37079 Goettingen, Germany. By use of this product the purchaser accepts the terms and conditions of all applicable Limited Use Label Licenses.

Trademark information

The owners of trademarks marked by "®" or "TM" are identified at http://www.iba-lifesciences.com/patents.html. Registered names, trademarks, etc. used in this $\ document, even \ when \ not \ specifically \ marked \ as \ such, \ are \ not \ to \ be \ considered \ unprotected \ by \ law.$