



Primary Mouse Whole Heart Dissociated Cells

Catalog Number: PC35136 **Components:** • Dissociated heart of E18 or P2 Mouse - >1 million viable cells

• 12 mls Culture Media- 12 ml of NbCardio: Cardiomyocyte growth media

Description: Fresh dissociated heart cells from E18 or P2 C57/Bl6 Mouse and media provided to initiate a culture of cardiomyocytes.

Shipping/Storage: Use dissociated cardiomyocytes immediately for highest cell yield; however, tissue can be stored for one week at 4-8°C. Dissociated heart is provided in 2 ml of Cardio Transport /Maintenance media.

Culture Protocol

Preparation (Room Temperature in a Sterile Hood)

1. Prepare substrate by coating slips or well plates with Gelatin (coating is 0.1% in filter sterilized DI water). Allow to sit for at least 3 hours to overnight at room temperature, wash once with DI water and allow to dry completely.
2. Fire polish the tip of a sterile 9" silanized pasteur pipette to an opening of ~0.5 mm
3. Aliquot 80 µl of Trypan Blue (Sigma: T8154) into a 0.5 ml tube for Step 6.

Cell Dispersal (Room Temperature in a Sterile Hood)

1. Allow 12 mL of NbCardio to come to room temperature.
2. Warm the 2 mL tube of dissociated cardiomyocytes in a 30°C water bath for 1 min.
3. With the silanized pasteur pipette, transfer the entire tube's contents to a sterile 15 ml centrifuge tube and triturate no more than 5 times. There should be no visible aggregation of cells in the media.
4. Spin 2000 rpm (~650 x G), 5 min. Discard supernatant leaving ~50 µl of Cardio Transport/ Maintenance media containing the pellet.
5. Disperse the pellet of cells (flick the bottom of the tube with a finger) and resuspend pellet in 1 ml NbCardio
6. Aliquot 20 µl of cell solution into the 0.5 ml tube containing 80 µl of Trypan Blue (1:5 dilution).
7. Count cells using a hemacytometer (calculate cells/ml).

Cell Plating (Room Temperature in a Sterile Hood)

1. Dilute cells with cardiac cell culture media (0.2 ml/cm²) and plate at 75,000 cells/cm² or desired concentration on 0.1% gelatin coated slips/flasks/wells.
2. Incubate 37°C, 5% CO₂, 9% O₂, 95% humidity (or ambient O₂)
3. After 24 hours, cardiomyocytes display spontaneous contracting and syncytial contracting at 120 hours
4. Change the medium with fresh, 37°C, CO₂ equilibrated cardiac cell media or media of choice every 2 days.
 - a. Around day 5 when the cells are confluent and beating in rhythm, we suggest switching from cardiomyocyte growth media to cardiomyocyte maintenance media

Viability Assay

1. Rinse cells twice with 37°C HBSS (0.2 ml/cm² of substrate).
2. Prepare dye mix from an acetone stock of 15 mg/ml fluorescein diacetate and an aqueous stock of 4.6 mg/ml propidium iodide, dilute 15 µl of each into 1.5 ml HBSS (1:100 dilution).
3. Add 20 µl of dye mix from step 2 to every 0.2 ml of HBSS added in step 1 (1:10 dilution).
4. After ~1 min count live cells using blue excitation appropriate for fluorescein fluorescence (green cells). Count dead cells with green excitation for propidium iodide fluorescence (small red nuclei).
5. Viability = (green cells/unit area)/(total cells plated/unit area) or Survival = green cells/(green + red cells)

FOR RESEARCH USE ONLY V5-06/2011

Neuromics' reagents are for in vitro and certain non-human in vivo experimental use only and not intended for use in any human clinical investigation, diagnosis, prognosis, or treatment. We disclaim all liability in connection with the use of the information contained herein or otherwise, and all such risks are assumed by the user.
